



ermes

CONNECT
YOUR PRODUCT
NOW

AREC Connect



JOINTLAB.COM
Jointlab S.r.l.

Via C. Treves 57, Trezzano sul Naviglio (Mi)
+39 0239310823 - info@frigolab.it

How to Connect to VELP Ermes Platform AREC Connect

In order to speed up the connection, be sure to get in contact with your IT specialist to get all your laboratory Wi-Fi connection details.

Requirements:


1. **Device:** PC, Tablet, Smartphone with Wi-Fi connection
2. **Browser:** Chrome, Safari, Firefox, Edge, UC browser

FIRST STEP | ACTIVATE YOUR VELP ERMES ACCOUNT

- Go to www.velp.com through your internet browser.
- In the Smart Lab menu click on "Configure Your VELP Ermes Account". Fill in the form and accept Terms and Conditions.
- You will receive an email containing your ID, Password and a link to proceed to VELP Ermes authentication at the email address that you entered during the registration.
- At the first login on VELP Ermes platform a new password will be requested: *remember to change it!*

SECOND STEP | CONNECT THE INSTRUMENT TO YOUR WI-FI NETWORK

- Turn on your AREC Connect
 - A. Press both knobs for 3 seconds; the display shows *TEMP LIMT*
 - B. Press the central button until *UIFI* appears; then press the right knob, rotate it to select *YES* and press it to confirm
 - C. Press the central button until *StAr AP* appears; then press the right knob, rotate it to select *YES* and press it to confirmThe Wi-Fi led starts blinking quickly

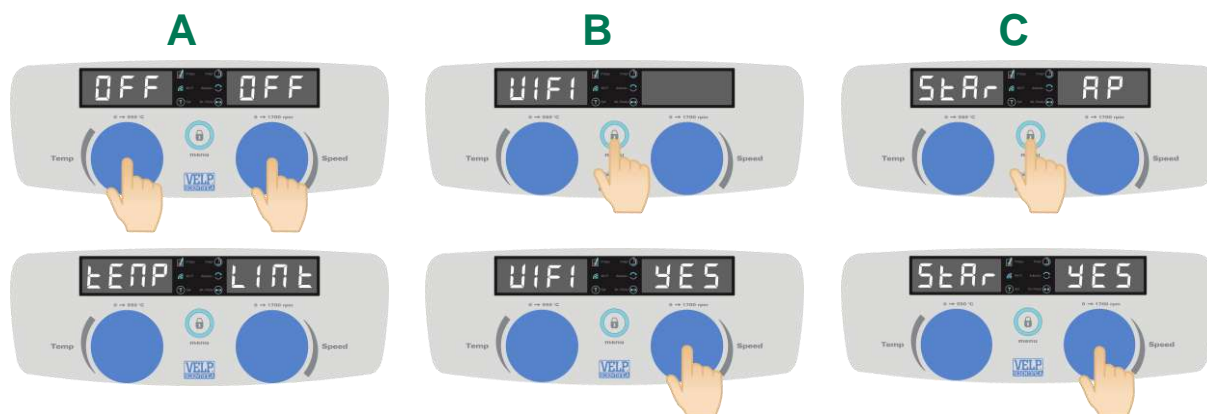

- From your device, select the VELP_AREC_(Product Code)_(Serial Number) network available on the Wi-Fi list.
- Open a web browser from your device (PC, tablet, smartphone), insert 192.168.4.1 into the address bar and press Enter. Use **admin** as user ID and **velp_ermes** as password.

Configure the laboratory Wi-Fi network parameters and press Push to confirm.

EXAMPLE:

Wi-Fi Name: (your Wi-Fi name)
Wi-Fi Password: (your Wi-Fi password)
Choose Auth: Wpa & WPA2 Personal
Choose Mode: Station

- Wait until the Wi-Fi led stops blinking.

NOTE: if the led is still blinking after 2 minutes, please try to repeat the procedure and check that the network parameters are correct



THIRD STEP | REGISTRATION ON VELP ERMES PLATFORM

A



Press both knobs for 3 seconds.

B




Press the central button until *OTC cFG* appears, then press the right knob.

C



On the display you will see "----".

D

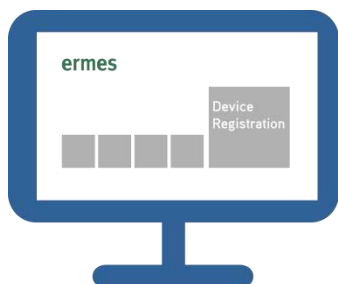

Login on VELP Ermes with your ID and Password and select "Device Registration".

E



Enter the AREC Connect product code, serial number (you can find this information on the silver label on the back of it) and purchase date. Press Next.

F




On the display of the AREC Connect will appear a four digits One Time Code (OTC). Enter this number, press Next and wait for connection.

G



Give a name to your AREC Connect and enter the address where your instrument is located. Press Next

H

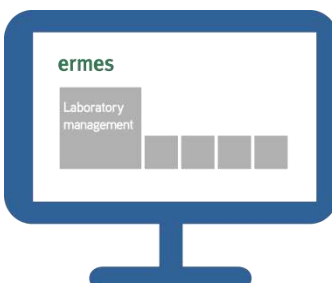

Add your AREC Connect to a laboratory or create a new one selecting "Laboratory Management".

I



Now from "Instruments" you can monitor and control your connected AREC Connect.



Always **Connected** and **Easy to Use**

Monitor and manage multiple instruments 24/7 from your device. Real time visualization of your analysis and of the instrument working conditions. Immediate notifications and alerts.



Safely Manage **Your Data**

VELP Ermes collects and stores your data with the maximum level of encryption following the highest cyber-security standards. You are now able to have access to your data anytime, anywhere. VELP Ermes will allow you to easily generate and share reports of your analyses.



Your Instrument **Always Updated**

Get software updates (OTA) immediately and benefit of all new features coming from VELP continuous product innovation.



Enhanced **Service Support**

Faster and smarter Service Support thanks to the direct contact with VELP specialists. With VELP Ermes we can guide you to get the maximum from your instrument.